

Why is Reversed-phase Flash Chromatography Use Increasing?

John R. Bickler, MA
Biotage, 10430 Harris Oaks Blvd., Suite C, Charlotte North Carolina 28269, USA

Introduction

Flash chromatography is the primary tool used by medicinal chemists for reaction mixture purification. Over the past 10 or so years, the use of reversed-phase flash chromatography for reaction mixture purification has increased dramatically due to both chromatographic and non-chromatographic reasons.

Some reasons for this switch from normal-phase include...

- » Synthesis of increasingly polar compounds
- » Synthesis of increasingly complex reaction mixtures
- » Conservation/environmental concerns/green chemistry
- » Cost savings
- » RxN mix analysis performed by HPLC-MS
- » Column reusability

In this poster, we will show an example of why chemists are increasingly using reversed-phase flash chromatography for purification of a complex reaction mixture.

Experimental Protocol

Reagents and Materials

Solvents used in study included: house DI water, hexane, ethyl acetate, acetonitrile and methanol (Reagents, Inc., Charlotte, NC).

Reagents used isatoic anhydride, benzyl amine, benzaldehyde, and acetic acid, all from Aldrich Chemical, Milwaukee, WI.

A Biotage[®] Initiator+ performed the synthesis.

A Biotage TLC plate (5x10 cm) was used for method development.

Flash chromatography was performed using a Biotage[®] Dalton 2000 system with an APCI ion source.

Normal-phase flash chromatography used a 5-gram, Biotage[®] Sfär HC column and Samplet[®] cartridge for dry loading.

Reversed-phase flash chromatography used a 6-gram Biotage[®] Sfär C18 flash column.

A Biotage[®] V-10 Touch dried the reaction mixture.

Organic Synthesis

Isatoic anhydride, benzyl amine, benzaldehyde, and acetic acid were mixed with acetonitrile and reacted to create 3-benzyl-2-phenyl-2,3-dihydro-4(1H)-quinazolinone, Figure 1. The reaction mixture concentration was 247 mg/mL.

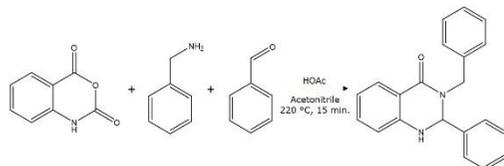


Figure 1. Microwave-assisted organic synthesis reaction in acetonitrile at 220 °C.

Thin-layer Chromatography

After synthesis, the reaction mixture was evaluated by TLC in 40% EtOAc in hexanes and separated a leading and trailing impurity away from the target product, Figure 2.



Figure 2. Reaction mixture TLC using 40% EtOAc/hexanes shows the product (blue) with a leading and trailing impurity.

Mass Analysis

The reaction mixture's detectable masses were determined using the Isolera Dalton 2000 with the solids analysis probe. Three m/z were found, 314.8 (product), 226.8, and 235.4, Figure 3.



Figure 3. Positive ionization APCI mass analysis finds three masses, 314.8 (product), 226.8 and 235.4

Flash Chromatography

A normal-phase linear gradient based on the TLC data (10-80% ethyl acetate) provided a partial separation as expected by the TLC results.

The flash method separated the lead impurity (m/z 235.4) from the other compounds but generated little resolution between the product and the trailing impurity (m/z 226.8), Figure 4.

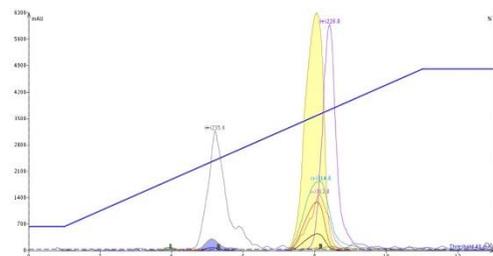


Figure 4. Reaction mixture normal-phase purification with mass identification. One impurity (m/z 235.4) is fully separated from the product but the other impurity (m/z 226.7) is not.

Reversed-phase

With normal-phase not providing an adequate purification result, I evaluated reversed-phase flash chromatography with a 50-80% acetonitrile in water gradient. The rationale being that reversed-phase's different separation mechanism

(partitioning vs. adsorption/desorption) may provide an improved purification result.

Reversed-phase flash chromatography improved separation with the m/z 226.8 impurity now eluting earlier than the product and the other impurity (m/z 235.4) eluting last, Figure 5. In fact, reversed-phase flash even resolved a previously un-detected impurity with a m/z 191.

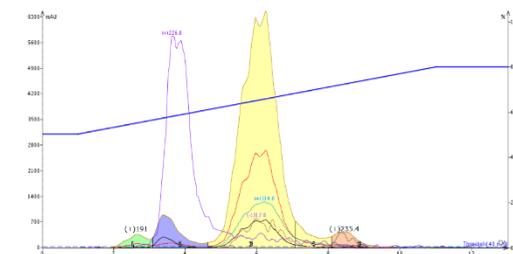


Figure 5. Reaction mixture reversed-phase flash purification separates both major impurities away from product. One impurity (m/z 226.8) elutes earlier than the product while the other impurity (m/z 235.4) elutes well after with this solvent system. Another impurity with m/z 191 is also separated.

Conclusions

For reaction mixtures not well-purified using normal-phase methods, reversed-phase flash chromatography often provides better separations because of differing separation mechanisms. This is a major reason why reversed-phase flash chromatography use is increasing in medicinal chemistry research labs.

Besides improved purification results, reversed-phase is greener and, generally, less expensive than normal-phase.