



## SPE-DEX 3000 Oil and Grease Spare Parts

To keep the extractor in good working conditions, the following items should be checked and replaced per the recommended frequency below.

Ref	Part #	Quantity for extractor	Description	Frequency
1	165-0712	1	Hydrophilic Solution, 250 mL	Perform every 60—90 Days to prevent Vapor Lock
2	48-2772	1 per Solvent Bottle	Service Kit, Solvent Bottle Filter	Replace every 12 months
3	50-0471	3 per System	Water Sample Inlet Valve	Inspect every 6 months, replace if worn or damaged
4	50-0667	1	Service Kit 1/3000 Solvent Rod Replacement Kit - 4 per Kit	Inspect once a month, replace if worn or damaged
5	02-0562	1	Black Rubber Stopper Plug for the 20 Liter Water-to-Waste Carboy	Inspect every 6 months, replace if worn or damaged
6	27-0042	5 for a 3000 System	Bottle - Plastic-Coated, 2.5 L (Solvent Delivery and Solvent-to-Waste)	Inspect Daily, replace if worn or damaged; replace
7	02-0006	1 per Solvent Bottle	Solvent Bottle Cap Gasket	Inspect every 6 months, replace if worn or damaged
8	50-0486-11	4	Solvent Delivery Lines from Bottles to Extractor	Inspect every 6 months, replace if worn or damaged
9	02-0216	1	Teflon® Gasket for the VOA Vial Adapter	Inspect every 6 months, replace if worn or damaged
10	50-0043-03	1	Solvent-to-Waste Manifold - (Top of Solvent-to-Waste Bottle)	Inspect every 6 months, replace if worn or damaged
11	50-0051-03	1	Water-to-Waste Manifold (Top of Water-to-Waste Carboy)	Inspect every 6 months, replace if worn or damaged
12	50-0485	1	Water-to-Waste and Solvent-to-Waste Lines Kit	Inspect every 6 months, replace if worn or damaged
13	50-0484	1	Gas Line, 3000 Series (Nitrogen Line)	Inspect every 6 months, replace if worn or damaged
14	50-0808-01	3 per System	SPE Disk Holder - 47 mm	Inspect every 6 months, replace if worn or damaged
15	50-0810-01	3 per System	SPE Disk Holder - 90 mm	Inspect every 6 months, replace if worn or damaged
16	30-0452-01	3 per System	"O" Ring on the bottom of Disk Holders	Inspect every 6 months, replace if worn or damaged



## SPE-DEX 3000 Oil and Grease Spare Parts

Continued

Ref	Part #	Quantity for extractor	Description	Frequency
17	31-0661	3 per System	Support screen for the 47-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
18	31-0662	3 per System	Support screen for the 90-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
19	31-0663	3 per System	Fine Mesh Screen for the 47-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
20	31-0665	3 per System	Fine Mesh Screen for the 90-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
21	02-0392	3 per System	Collection Adapter	Replace Yearly
22	30-0441	3 per System	"O" Ring used on the Collection Adapter	Replace Yearly
23	48-2605	3 per System	Sample-Collect Check Valve Kit	Inspect every 6 months, replace if worn or damaged
24	27-0834-01	3 per System	Separatory Funnel - 125-mL, 19/22 Taper	Inspect every 6 months, replace if worn or damaged
25	22-0687	3 per System	Blue Clamp used to secure the collection vessel to the Collection Adapter	Inspect every 6 months, replace if worn or damaged
26	02-1984	1	Check Valve Release Tool	Inspect every 6 months, replace if worn or damaged